*USB Microscope M3*

# Hardware Manual

English/Japanese

**SCALAR**  
IMPRESSIVE WORKS

## Operating environment

### Required environment

Preinstalled Windows XP SP2 or later  
Windows Vista

USB Port : Factory installed USB2.0 or USB1.1 Port

CPU : Pentium 4 or equivalent

### Cautions


- If two or more USB devices are connected to a single computer, the product may not function properly, depending on the type of USB device being used simultaneously.
- Product compatibility is not guaranteed if this product is used through a hub.
- \* Product compatibility is not guaranteed for all computers, even having the recommended operating environments.

\* Windows® is registered trademark of Microsoft Corporation.

\* Other Product names are registered trademarks and trademarks of their respective holders.

## USB Microscope Schematic

### Parts of the USB Microscope

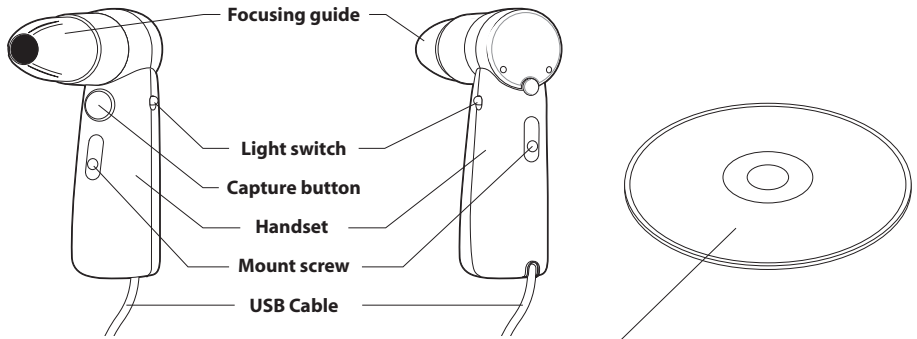**CD-ROM (Application Software for Windows®)**

\* You can download the latest version at [www.scalar.co.jp/](http://www.scalar.co.jp/)

### Specifications

Imager	1/3 inch high resolution 2Mp color sensor 1,920,000 effective pixels	
Interface	USB 2.0 or USB 1.1. USB 2.0 required for high resolution mode SXGA (1280x1024), and higher frame rates.	
Magnification	Varies based on attached lens.	
Weight	About 200g with lens unit attached.	
Power consumption	About 400mA, when using a 5V light source	
Length of Cable	About 1.8m (6 ft)	
Capture Button	Initiates still, movie, and time lapse capture in Scalar software.	
Light Switch	3-position on/off switch for lens LEDs, A-OFF-B	
Camera Mount	standard tripod mount, 1/4" - 20 threads	
Lens Release Button	Disengages lens lock pin for lens removal.	
Video Preview	<b>USB 2.0 High Speed</b> 30 fps @ 320x240 pixels (QVGA) 15 fps @ 640x480 pixels (VGA) 3.75 fps @ 1280x1024 pixels (SXGA)	<b>USB 1.1 Full Speed</b> ~4fps@320x240pixels (QVGA) ~1fps@640x480pixels (VGA) High resolution mode (SXGA) not available.
Video Controls	Brightness, Contrast, Saturation, and Sharpness. Light Source/White Balance options include M3 LED, Incandescent, Fluorescent, Daylight, and Auto.	
Still Capture	320x240 pixels (QVGA) 640x480 pixels (VGA) 1280x1024 pixels (SXGA) A USB 1.1 port may be used but only 640x480(VGA) and 320x240 (QVGA) modes are supported and at a much reduced frame rate.	
Vodeo Capture	Same as the video preview.	
Driver	USB Video Class, USB Human Interface Device (HID) Class. WindowsXP SP2 or later, WindowsVista	
System Requirements	Windows XP Service Pack 2 or later. Factory installed USB 2.0 or USB 1.1 port. USB 2.0 port recommended. A USB 1.1 port may be used but only 640x480(VGA) and 320x240 (QVGA) modes are supported and at a much reduced frame rate. Pentium 4 or equivalent	
Operating Temperature Range	5°C~40°C	
Storage Temperature Range	-10°C~60°C (without condensation)	

## Software Installation

Install application software from attached CD-ROM into your PC.  
Refer to Software Manual in CD-ROM for operation instruction.

## Connecting the USB Microscope

Note: The USB Microscope will work only with computers that support the USB interface.

Once your computer has finished starting up, connect the USB Microscope cable to a USB port on your computer.

**Note:** For the location of the USB port(s) on your computer, refer to your computer's user's manual.

To ensure that the USB Microscope functions normally, do not connect or disconnect the USB cable while the system is starting or closing.

To avoid unstable system software functioning, do not connect or disconnect a USB cable while an application is opening or closing.

When disconnecting the USB Microscope with your computer turned on, observe the steps 1 and 2 following. Disconnecting improperly may cause the computer to operate abnormally.

Step 1 is not necessary if the computer is turned off. Note that the computer may not recognize the microscope if the USB cable is disconnected and then reconnected while the computer is in power-saving mode.

The correct way to disconnect the microscope:

1. Close any software currently using the USB Microscope.
2. Disconnect the USB cable.



Never pull on the cable portion to disconnect the USB cable. Always grasp the connector itself when disconnecting.

See "Software Manual" in the CD-ROM for using the application program.

## Using the USB Microscope

### Focusing guide


Select either of the following two positions for the focusing guide of the M50 lens unit or M30N lens unit supplied with the product: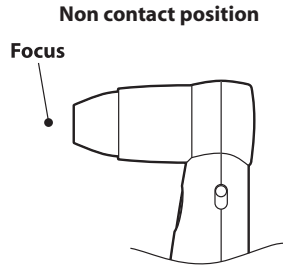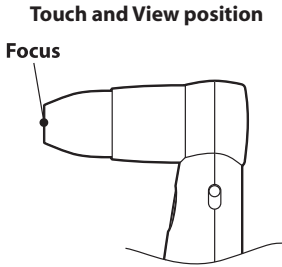

**Touch and View Position**-The tip of the focusing guide is placed in direct contact with the object.

**Noncontact Position**-This position is selected to observe objects with significant surface roughness, or when the USB Microscope is attached to a stand for observations.

Observations are made without placing the focusing guide into contact with the object.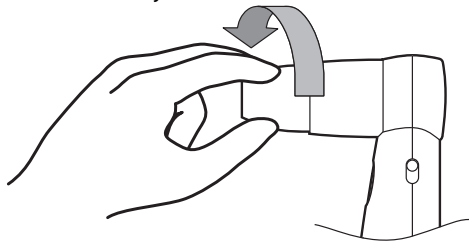

Set to the Touch and View position by turning the lens unit counterclockwise until you hear a click.

### Lighting switch

Slide the switch to turn the light on.


A: Reflection mode (for every lens unit)

Off: Light off


B: Non Reflection mode (only for M30N lens unit)

\* Non Reflection mode only works with M30N Lens unit.

\* Please turn off the light when you do not use the microscope for a long time.

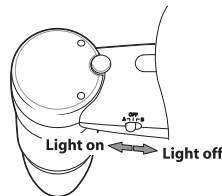
## Changing the Lens-unit
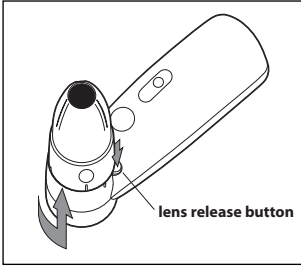
### 1. Detaching the Lens-unit

Turn the lens base counterclockwise while pressing the Lens Release Button until the hash mark on the Main Unit body and a dot on the lens line up.

**NEVER TRY TO REMOVE A LENS WITHOUT USING THE LENS RELEASE BUTTON !**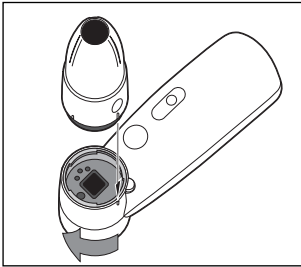

### 2. Attaching the Lens-unit

Line up the hash mark on the Main Unit body and dot on the lens, insert the lens unit, and turn clockwise until the lens base is tight.

## Handling Notes

### ■ Avoid positioning the USB Microscope at the following locations

To avoid damage to the USB Microscope, avoid placing it in any of the following locations, either for use or for storage.

Locations subject to unusually high temperatures

Automobile interiors during the summertime, especially with windows closed. Such heat may deform or damage the USB microscope.

Locations in direct sunlight or near heating devices

The heat found in such environments may deform or damage the USB Microscope.

Locations subject to sandy particles, including sandy soil or beaches

Near beaches, on sandy soil, or in any place subject to sand-like dust, take adequate measures to protect the USB Microscope. Without protection, such environments may damage the USB Microscope beyond repair.

Do not keep the USB Microscope positioned with its lens directed toward the sun.

The lens may be damaged by scorching. Exercise caution when placing the USB Microscope close to a window or outdoors.

### ■ Operating Notes

Do not disconnect the USB cable of the USB Microscope while an application is running. Doing so can lead to system errors or instability.


Do not place any object on the USB Microscope.

Avoid sudden changes in ambient temperature. Do not use the USB Microscope immediately after moving it from a cold place to a warm place, or after sudden increase in room temperature. Such changes may produce condensation inside the USB Microscope. Allow the unit to stand for a while before use.

### ■ Transportation

Retain the original packing in case you need to transport the USB Microscope in the future.

### ■ Handling of the CMOS image sensor section

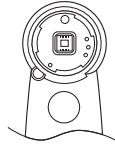The CMOS image sensor section is exposed when you remove the lens-unit from the handset.

Never touch the exposed CMOS image sensor section.

### ■ Handling of the CD-ROM

To protect data and information recorded on the CD-ROM, observe the following precautions.

Avoid scratching the CD-ROM surface.

When handling the CD-ROM, avoid touching the blank side. Data is recorded on this side.

Avoid storing the CD-ROM in dusty locations or in direct sunlight, near heat sources, or in extremely humid locations.

Take care to avoid spilling liquid on the CD-ROM surface.

Store the CD-ROM in a suitable storage case.


### ■ In case of problems or malfunction

Please consult with your dealer or distributor if the USB Microscope malfunctions, or if its interior becomes contaminated with foreign particles.

## Maintenance


Clean the exterior of the USB Microscope with a soft, dry cloth. In cases of heavy soiling, wipe the surface with cloth moistened with a dilute, neutral detergent. Never use thinner, benzene, or alcohol - such materials will damage the surface finish.

### Dust on the CMOS image sensor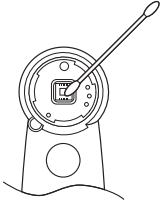


Dust on the CMOS image sensor section may be removed via the aperture with a blower or dust gun, or carefully removed with a suitable applicator.

### Dust on the lens-unit

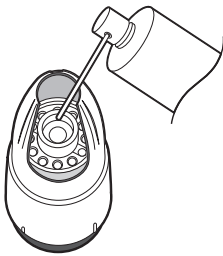Dust on the lens-unit surface may be removed via the aperture with a dust gun.

### Dirt on the CD-ROM

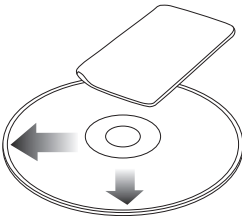If its surface is fouled by finger prints or dust, the CD-ROM may not be read reliably. Keep the surface clean at all times.

- Using a soft cloth, lightly wipe the CD-ROM surface from the center to the outer periphery.
- Clean heavily-soiled CD-ROM surfaces by wiping with a water-moistened cloth. Wipe dry with a dry cloth.
- Do not use benzene, record disk cleaner, or antistatic agents - they may damage the CD-ROM.

## Troubleshooting Guide - Quick solutions to save your time

### ■ The USB Microscope does not work. (The Power indicator does not lighten.)

- Check to make sure that the USB cable is plugged into the USB port. (See P3)

### ■ The application program is not on.

- Check to make sure that the USB cable is plugged into the USB port. (See P3)

### ■ The application program is on, but there is no picture.

- Check to make sure that the lamp of the lens-unit lightens rightly. (See P4)

### ■ The display screen shows incorrect colors.

- Adjust the appropriate items in Video settings in the application program. (See Software Manual)

### ■ Captured still images and animation images are indistinct on the display screen.

- Images captured under fluorescent lighting may appear on screen together with lamp-light reflection components.
- Captured images containing dark spots or areas may result in red or green color noise.
- The CMOS image sensor section may be fouled. -Clean the CMOS image sensor section. (See P6)
- The lens-unit is fouled. -Clean the lens-unit. (See P6)
- The image is out of focus. -Set the Focusing guide rightly. (See P4)

### ■ Frame skipping errors occur when capturing video.

- Running more than one application at a time may exceed your computer's processing capabilities. Close all other software.
- The computer's power-management function has been activated. Check the speed setting for the CPU.

For any inquiry, please contact your nearest Scalar reseller or [sales@scalar.co.jp](mailto:sales@scalar.co.jp)  
[www.scalar.co.jp/english](http://www.scalar.co.jp/english)

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**Scalar Corporation**